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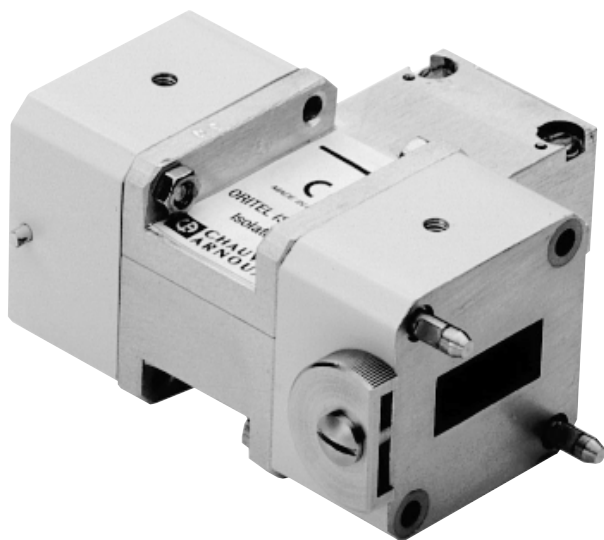
Meaning of the symbol

CAUTION! Consult the operating manual before using the component.

In this operating manual, the instructions preceded by this symbol may cause personal injury or damage the instrument and installations if they are not scrupulously abided by or carried out.

You have just bought an **ORITEL ISO 100 ferrite isolator** and we thank you for the interest you show for our products.

In order to get the best use out of your component:

- carefully **read** this operating manual,
- **follow** the precautions for use.

 PRECAUTIONS FOR USE

- Take care that the isolator is fitted in the correct direction, with the arrow pointing towards the load.
- All instruments hooked up to the bench and connected to the AC network shall be earthed.
- No voltage of more than 30 V in relation to the earth shall be present on the bench.

TO ORDER

ORITEL ISO 100 ferrite isolator P01.2753.08

Delivered with this operating manual.

Replacement part

ORITEL AFR 100 EASYFIX quick fastening lug P01.2753.01

CONTENTS

1. PRESENTATION	7
2. DESCRIPTION	7
3. USE	8
3.1 Summary of the components making up the ORITEL BDH R100 basic educational bench ...	8
3.2 How to use the EASYFIX quick fastening lug	8
4. CHARACTERISTICS	9
5. MAINTENANCE	9
5.1 Cleaning	9
5.2 Metrological inspection	9
6. GUARANTEE	9
7. ANNEX	22
7.1 General view of the ORITEL ISO 100	22
7.2 Assembling the EASYFIX quick fastening lug adaptor on the ORITEL ISO 100 isolator flanges	22
7.3 Using the EASYFIX quick fastening lug	23

1. PRESENTATION

The **ORITEL ISO 100 ferrite isolator** is a microwave component specially designed for the ORITEL BDH R100 educational bench, manufactured using a rectangular waveguide in accordance with the R100/WR90 standard (covering the frequency range between 8.5 and 9.6 GHz).

It enables the OSG 100 Gunn oscillator to be protected against variations in load impedance and against cases of mismatching caused by the MOD 100 pin diode modulator, with the effect of these mismatches being to cause frequency and power variations in the Gunn source.

The isolator has the distinctive feature of letting energy through with only a small insertion loss in the oscillator → load direction and brings about a high degree of attenuation in the opposite direction.

2. DESCRIPTION

See § 7.1

- ① Arrow indicating the direction in which the isolator should be fitted.
- ② Knurled tightening knob for the *EASYFIX* quick fastening lug.

3. USE

The ferrite isolator, used in all the manipulations described in the operating manual relevant to the ORITEL BDH R100 educational bench, **shall be correctly oriented, with the arrow indicating the load**. The components below are generally included in the standard supplies delivered with the ORITEL BDH R100 educational bench.

To use these components, refer to their respective operating manuals.

3.1 Summary of the components making up the ORITEL BDH R100 basic educational bench

ORITEL OSG 100 Gunn diode oscillator	P01.2753.07
ORITEL ISO 100 ferrite isolator	P01.2753.08
ORITEL MOD 100 pin diode modulator	P01.2753.09
ORITEL ATM 100 micrometer-adjustable variable attenuator	P01.2753.10
ORITEL OND 100 curve wavemeter	P01.2753.11
ORITEL LAF 100 measuring line	P01.2753.12
ORITEL ADZ 100/3 three-screw impedance adaptor	P01.2753.13
ORITEL TGN 100 waveguide-coaxial transition element	P01.2753.14
ORITEL DEN 100 coaxial detector	P01.2753.15
ORITEL CHG 100 adapted load	P01.2753.16
ORITEL CC 100 short circuit plate	P01.2753.17
ORITEL SUP 100 waveguide support	P01.2753.18

■ Other components (R100/WR90 standard waveguide)

ORITEL AFR 100 «EASYFIX» quick fastening lug adaptor	P01.2753.01
ORITEL RS 100 bench support rail	P01.2753.03
ORITEL MH 600 microwave milliwattmeter	P01.2501.01
ORITEL ST 613 thermocouple probe (for ORITEL MH 600)	P01.2851.01
ORITEL CF 204A power supply unit for Gunn diode oscillator	Please consult us
ORITEL IR 205 SWR indicator	P01.2705.01
ORITEL RD 100 displacement sensor for measuring line	P01.2753.02
ORITEL ANC 100/15 (15 dB) horn antenna	P01.2753.04
ORITEL CGX 100/20 waveguide cross coupler	P01.2753.05
ORITEL IRIS 20-30 20 and 30 dB) coupling irises	P01.2753.06

To use these components, refer to their respective operating manual.

3.2 How to use the EASYFIX quick fastening lug

See diagram § 7.3

- Arrange the two elements to be assembled, taking care to correctly place the centring studs ① opposite the holes ②
- Simultaneously tighten the two knurled knobs ③

4. CHARACTERISTICS

Frequency: 8.5 to 9.6 GHz

Insertion loss: < 1 dB

Isolation: > 20 dB

SWR: ≤ 1.25

Waveguide: R100

Flange: UBR 100

Weight: 650 g (with 2 *EASYFIX* quick fastening lug adaptors)

Dimensions: 99 x 46 x 87 mm

Electromagnetic compatibility:

Compliance with NF EN 61326-1 standard, 1998 edition.

5. MAINTENANCE

 **For maintenance, use only specified spare parts. The manufacturer will not be held responsible for any accident occurring following a repair done other than by its After Sales Service or approved repairers.**

5.1 Cleaning

Before assembling the educational bench's microwave components, make sure that no dust is present inside the waveguide.

CAUTION!

Do not use a pressurised air jet for fear of damaging certain components.

Do not use solvent, and take precautions when dusting.

5.2 Metrological inspection

 **It is essential that all measuring instruments are regularly calibrated.**

For checking and calibration of your instrument, please contact our accredited laboratories (list on request) or the Chauvin Arnoux subsidiary or Agent in your country.

■ **Repairs under or out of guarantee**

Please return the product to your distributor.

6. GUARANTEE

Our guarantee is applicable for **twelve months** after the date on which the equipment is made available (extract from our General Conditions of Sale, available on request).

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Deutschland : CA GmbH - Straßburger Str. 34 - 77694 Kehl / Rhein - Tel : (07851) 99 26-0 - Fax : (07851) 99 26-60
Schweiz : CA AG - Einsiedlerstrasse 535 - 8810 Horgen - Tel : (01) 727 75 55 - Fax : (01) 727 75 56
UK : CA UK Ltd - Waldeck House - Waldeck road - Maidenhead SL6 8br - Tel : (01628) 788 888 - Fax : (01628) 628 099
USA : CA Inc - 99 Chauncy Street - Boston MA 02111 - Tel : (617) 451 0227 - Fax : (617) 423 2952
USA : CA Inc - 15 Faraday Drive - Dover NH 03820 - Tel : (603) 749 6434 - Fax : (603) 742 2346